



*Suggestions or questions to:
Rebecca Williams, Director
Cornell Imaging, BRC
rw36@cornell.edu*

Antifade mounting medium recipe

Ingredients:

10X PBS stock solution

Glycerol

Dimethyl sulfoxide (DMSO)

n-propyl gallate (NPG, Sigma P3130)

Directions:

- 1) Prepare a stock solution of 20%(w/v) NPG in DMSO.
- 2) Mix 1 part of 10X PBS with 9 parts of glycerol.
- 3) Slowly add 0.1 part 20% NPG stock solution dropwise with rapid stirring.