

Metamorph Installation Instructions (For Windows only)

(Note: if you install this program in the Program Files location rather than the default location, you will need to run it in Administrator mode under Windows 7.)

- 1) Go to the Imaging FileShare.
- 2) In VIZ\Metamorph\METAMORPH_FOR_OLYMPUS, go either to the Setup or the Setup-x64 directory, depending on whether your computer has a 32- or 64-bit operating system. (If you don't know, use the 64-bit option. It will not install if inappropriate.)
- 3) Double click on setup.exe
- 4) Click Next and Accept the license terms.
- 5) Do not Enable Sentinel Component Installation and keep on clicking Next.
- 6) At the Hardware Key Confirmation menu, choose Network Key Server and in the Server Address field type:

keyserver2.biotech.cornell.edu
- 7) Click the Read System ID from Key button. (System ID 25228 should show up.)
- 8) Click Next.
- 9) At the User Information and Codes menu: For Retrieve Codes, click Using the Internet. Lots of letters and numbers should appear. Click Next.
- 10) Click Install. Click OK. Click Finish. (You may need to restart Windows.)
- 11) You should be done. Open Metamorph Basic Offline.